

Selectin Biosciences Inc.

Endoglycosidase H

Endoglycosidase H [Endo- β -N-acetylglucosaminidase H, EC 3.2.1.96] cleaves asparagine-linked oligomannose and hybrid, but not complex, oligosaccharides from glycoproteins (see Figure). It cleaves between the two N-acetylglucosamine residues in the diacetylchitobiose core of the oligosaccharide, generating a truncated sugar molecule with one N-acetylglucosamine residue remaining on the asparagine. In contrast, PNGase F removes the oligosaccharide intact. Detergent and heat denaturation may increase the rate of cleavage for some glycoproteins.

Endoglycosidase H is produced from a *Streptomyces plicatus* clone.

Product Code: GE 44

Specifications

Activity: 40 U/mg, ~5 U/ml

Storage: Store at 4°C. Do Not Freeze.

Formulation: The enzyme is provided as a sterile solution in 20 mM Tris HCl pH 7.5, 50 mM NaCl 1 mM EDTA.

Stability: Stable at least 12 months when stored properly. Several days exposure to ambient temperatures will not reduce activity.

Product Description

Molecular weight: 29,000 daltons

Purity: Endoglycosidase H is tested for contaminating protease as follows; 10 μ g of denatured BSA is incubated for 24 hours at 37°C with 2 μ l of enzyme. SDS-PAGE analysis of the

treated BSA shows no evidence of degradation. The production host strain has been extensively tested and does not produce any detectable glycosidases.

Specificity: Asparagine-linked hybrid or high mannose oligosaccharides.

Assay

One unit of Endo H activity is defined as the amount of enzyme required to catalyze the release of N-linked oligosaccharides from 1 μ mole of denatured Ribonuclease B in 1 minute at 37°C, pH 5.5. Cleavage is monitored by SDS-PAGE (cleaved Ribonuclease B migrates faster).

Reagents

- 5X PNGase buffer 5.5- 250 mM sodium phosphate pH 5.5.
- Denaturation solution- 2% w/v sodium lauryl sulfate, 1 M β -mercaptoethanol.

Suggestions for Use

Procedure for deglycosylation

1. Add up to 200 μ g of glycoprotein to Eppendorf tube. Adjust to 37.5 μ l final volume with deionized water.

2. Add 10 μ l 5X Endoglycosidase H Buffer and 2.5 μ l of Denaturation Solution (SDS/-ME). Heat at 100°C for 5 minutes.

NOTE: It is not necessary to add Triton X-100. SDS will not inactivate Endoglycosidase H.

Selectin Biosciences Inc.

3. Cool. Add 2.0 µl of Endoglycosidase H to the reaction. Incubate 3 hours at 37°C.

If SDS or heat denaturation is omitted, it may be necessary to increase incubation time.

References

1. Robbins P. W., D. F. Wirth and C. J. Hering. Expression of the Streptomyces enzyme endoglycosidase H in Escherichia coli. **Biol Chem 256**:10640-10644 (1981).
2. Robbins P. W., R. B. Trimble, D. F. Wirth, C. Hering, F. Maley, G. F. Maley, R. Das, B. W. Gibson, N. Royal and K. Biemann. Primary structure of the Streptomyces enzyme endo-beta-N-acetylglucosaminidase H. **J Biol Chem 259**:7577-7583 (1984).
3. Trimble R. B., A. L. Tarentino, G. E. Aumick and F. Maley. Endo-beta-N-acetylglucosaminidase L from Streptomyces plicatus. **Methods Enzymol 83**:603-610 (1982).
4. Trimble R. B. and F. Maley. Optimizing hydrolysis of N-linked high-mannose oligosaccharides by endo-beta-N-acetylglucosaminidase H. **Anal Biochem 141**:515-522 (1984).
5. Trimble R. B., R. J. Trumbly and F. Maley. Endo-beta-N-acetylglucosaminidase H from Streptomyces plicatus. **Methods Enzymol 138**:763-770 (1987).
6. Trumbly R. J., P. W. Robbins, M. Belfort, F. D. Ziegler, F. Maley and R. B. Trimble. Amplified expression of Streptomyces endo-beta-N-

acetylglucosaminidase H in Escherichia coli and characterization of the enzyme product. **J Biol Chem 260**:5683-5690 (1985).

Selectin Biosciences Inc warrants that the above product conforms to the specifications described herein. Should the product fail for reasons other than through misuse Selectin Biosciences Inc. will, at its option, replace free of charge or refund the purchase price. This warranty is exclusive and Selectin Biosciences Inc. makes no other warrants, expressed or implied, including any implied conditions or warranties of merchantability or fitness for any particular purpose. Selectin Biosciences Inc. shall not be liable for any incidental, consequential or contingent damages.

This product is intended for in vitro research only.

REVISION 6/15/15