

Selectin Biosciences Inc.

Endoglycosidase F1

Endoglycosidase F1 [Endo- β -N-acetylglucosaminidase F1, EC 3.2.1.96] cleaves asparagine-linked or free oligomannose and hybrid, but not complex, oligosaccharides (see figure 1). Core fucosylation reduces the activity by 50 fold. Endoglycosidase F1 will hydrolyze sulfate containing high-mannose chains. It cleaves between the two N-acetylglucosamine residues in the diacetylchitobiose core of the oligosaccharide, generating a truncated sugar molecule with one N-acetylglucosamine residue remaining on the asparagine. In contrast, PNGase F removes the oligosaccharide intact.

Endoglycosidase F1 is less sensitive to protein conformation than PNGase F and is therefore more suitable for deglycosylation of native proteins. However for optimal results, denaturation of the glycoprotein is recommended.

Endoglycosidase F1 is isolated from a strain of *E. coli* expressing a cloned gene from *Elizabethkingia miricola*.

Product Code: GE 47

Specifications

Activity: ≥ 16 U/mg, ≥ 17 U/mL

Storage: Store at 4°C. Do not freeze.

Formulation: The enzyme is provided as a sterile solution in 20 mM Tris HCl, pH 7.5.

Stability: Stable at least 12 months when stored properly. Several days exposure to ambient temperatures will not reduce activity.

Product Description

Molecular Weight: 32,000 daltons

Purity: Endoglycosidase F1 is tested for contaminating protease as follows; 10 μ g of denatured BSA is incubated for 24 hours at 37°C with 2 μ L of enzyme. SDS-PAGE analysis of the treated BSA shows no evidence of degradation.

The production host strain has been extensively tested and does not produce any detectable glycosidases.

Specificity: Asparagine-linked hybrid or free hybrid or high mannose oligosaccharides.

Assay

One unit of Endoglycosidase F1 activity is defined as the amount of enzyme required to catalyze the release of N-linked oligosaccharides from 1 μ mole of denatured Ribonuclease B in 1 minute at 37°C, pH 5.5. Cleavage is monitored by SDS-PAGE (cleaved Ribonuclease B migrates faster).

Reagents

- 5X Reaction buffer 5.5 – 250 mM sodium phosphate pH 5.5
- Denaturation Solution: (2% SDS, 1 M β -mercaptoethanol [β -ME])
- Triton X-100 solution, 15%

Suggestions for Use

Procedure for Deglycosylation

1. Add up to 200 μ g of glycoprotein to Eppendorf tube.

Selectin Biosciences Inc.

2. Add deionized water to a total of 33 μ L.
3. Add 10 μ L 5X Reaction Buffer, 5.5.
4. Add 2.5 μ L of Denaturation Solution. Heat at 90°C for 10 minutes.
5. Cool to room temperature and add 2.5 μ L Triton X-100 solution.
6. Add 2 μ L of Endoglycosidase F1. Incubate 1 hour or more at 37°C.
7. Monitor cleavage by SDS-PAGE.

For digestion of native proteins, add water to a total volume of 38 μ L and omit steps 4 and 5. Increase incubation time appropriately.

References

1. Maley P., R. B. Trimble, A. L. Tarentino and T. H. Plummer Jr. Characterization of glycoproteins and their associated oligosaccharides through the use of endoglycosidases. **Anal Biochem** **180**:195-204 (1989).
2. Plummer, T. H. Jr, A. W. Phelan and A. L. Tarentino. Porcine fibrinogen glycopeptides: substrates for detecting endo-beta-N-acetylglucosaminidases F2 and F3. **Anal Biochem** **235**: 98-101 (1996).
3. Reddy A., B. G. Grimwood, T. H. Plummer Jr and A. L. Tarentino. High-level expression of the Endo-beta-N-acetylglucosaminidase F2 gene in E.coli: one step purification to homogeneity. **Glycobiology** **8**:633-636 (1998).
4. Tarentino, A. L., C. M. Gomez and T. H. Plummer Jr. Deglycosylation of Asparagine-Linked Glycans by Peptide:N-Glycosidase F. **Biochemistry** **24**:4665-4671 (1985).
5. Tarentino A. L., G. Quinones, W. P. Schrader, L. M. Changchien and T. H. Plummer Jr. Multiple endoglycosidase (Endo) F activities expressed by *Flavobacterium meningosepticum*. Endo F1: molecular cloning, primary sequence, and structural relationship to Endo H. **J Biol Chem** **267**:3868-3872 (1992).
6. Tarentino A. L., G. Quinones, L. M. Changchien, and T. H. Plummer Jr. Multiple endoglycosidase F activities expressed by *Flavobacterium meningosepticum* endoglycosidases F2 and F3: Molecular cloning, primary sequence, and enzyme expression. **J Biol Chem** **268**(13):9702-9708 (1993).
7. Tarentino A. L. and T. H. Plummer Jr. Substrate specificity of *Flavobacterium meningosepticum*: Endo F2 and endo F3: purity is the name of the game. **Glycobiology** **4**:771-773 (1994).
8. Tarentino, A. L. and T. H. Plummer Jr. Enzymatic deglycosylation of asparagine-linked glycans: purification, properties and specificity of oligosaccharide-cleaving enzymes from *Flavobacterium meningosepticum*. **Methods in Enzymology** **230**:44-57 (1994).
9. Tarentino A. L., G. Quinones and T. H. Plummer Jr. Overexpression and purification of non-glycosylated recombinant endo-beta-N-acetylglucosaminidase F3. **Glycobiology** **5**:599-601 (1995).

Selectin Biosciences Inc.

10. Trimble, R. B. and A. L. Tarentino.
Identification of Distinct
Endoglycosidase (Endo) Activities in
Flavobacterium meningosepticum: Endo
F1, Endo F2 and Endo F3. **J. Biol Chem**
266:1646-1651 (1991).

Selectin Biosciences Inc warrants that the above product conforms to the specifications described herein. Should the product fail for reasons other than through misuse Selectin Biosciences Inc. will, at its option, replace free of charge or refund the purchase price. This warranty is exclusive and Selectin Biosciences Inc. makes no other warrants, expressed or implied, including any implied conditions or warranties of merchantability or fitness for any particular purpose. Selectin Biosciences Inc. shall not be liable for any incidental, consequential or contingent damages.

.

This product is intended for in vitro research only.

REVISION 7/31/12

Selectin Biosciences Inc.

Figure 1 - Cleavage of oligosaccharides by Endo F1 and PNGase F

Man - Mannose; GlcNAc - N-acetylglucosamine

